OCT 15 1984

**NOTICE OF APPROVAL**

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Company: Fisher Scientific  
Division of Allied Canada Inc.  
112 Colonnade Road  
NEPEAN, Ontario  
K2E 7L6

Société: Fisher Scientific  
Division of Allied Canada Inc.  
112 rue Colonnade  
NEPEAN, Ontario  
K2E 7L6

Manufacturer: Mettler Instrument AG  
C H-8606 Greifensee  
Switzerland

Fabricant: Mettler Instruments AG  
C H-8606 Greifensee  
Switzerland

Type of Device: Electronic top loading  
balance.

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NOTE: \*Delta Range - this is a mode that increases the increment by a factor of 10 and reduces capacity by the same factor.

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PE models with a Delta Display whereby the last decimal place is blanked out on initial application of weight but as the weight settles it re-appears. This function speeds up the weighing cycle.

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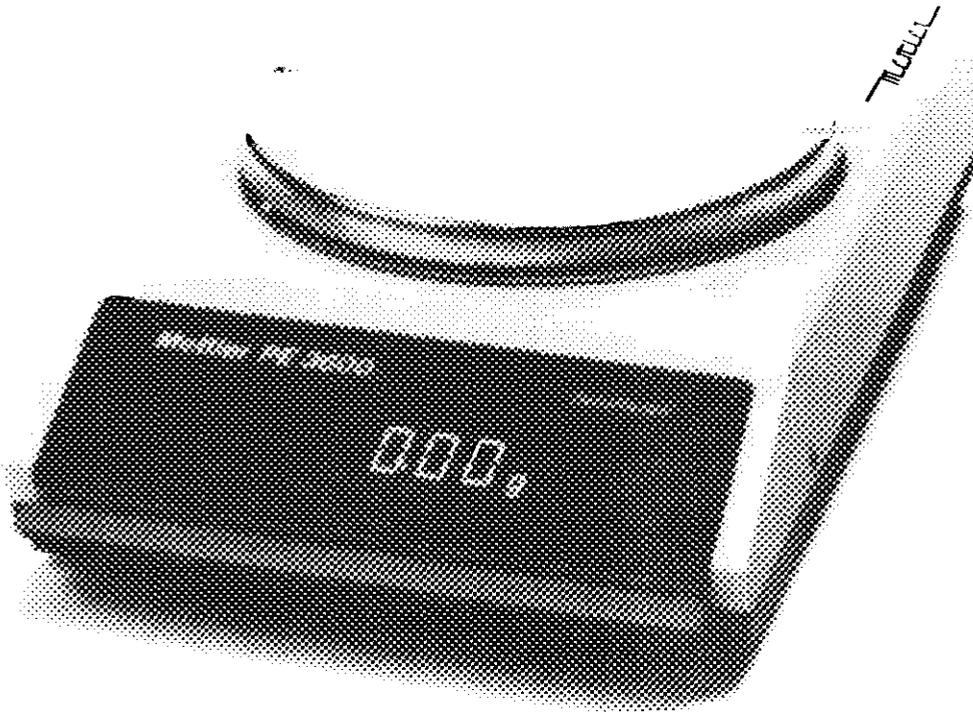
Accessories: Compatible accessories that meet the requirements and perform in accordance with the Weights and Measures Act, Regulations and Specifications may be used.

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.../2

- 2 -

Application: For exclusive use in weighing precious metals and other commodities of comparable value.

Description: These are top loading electronic balances. Weight is sensed using an electromagnetic force compensation principle. All models are fitted with lockable levelling feet and a bullseye level. These balances are fitted with + and - indication to the left of the weight display; when both are off, a zero display indicates that the zero position is accurate to  $\pm \frac{1}{4}$  of an increment.

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Description: Continued

There is only one operator control for these balances which controls, ON/OFF, zero/tare and calibrate.

To switch ON - press bar

To switch OFF - lift bar

To zero - press bar. It operates to full scale capacity.

To tare - press bar. It operates to full scale capacity.

Blanking for over capacity is indicated by a series of flat bars on the weight display.

These balances have one display only, they must be labelled with instructions that state "THIS BALANCE MUST BE LOCATED SO THAT THE WEIGHT DISPLAY IS VISIBLE TO ALL PARTIES OF ANY TRADE TRANSACTION" (or words to that effect).

These balances must be installed on a solid, stable surface free from vibration and in an environment free from air currents.

Zero exceeds 4% of capacity.

Tare is not identified nor is there visual confirmation when the tare mode is operating.

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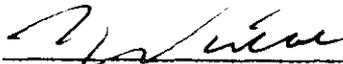
Conditions of Approval: Approval is granted under the Weights and Measures Act, S.C. 1970-71-72, chapter 36, and the Weights and Measures Regulations C.R.C.c. 1605 for use in Canada under the general conditions of the said Regulations, and under any special conditions listed above.

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  
W.R. Virtue

Chief  
Legal Metrology Laboratories

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