NOTICE OF APPROVAL

Issued by statutory authority of the Minister of Industry for:

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NOTE: This approval applies only to devices, the design, composition, construction and performance of which are, in every material respect, identical to that described in the material submitted, and that are typified by samples submitted by the applicant for evaluation for approval in accordance with sections 14 and 15 of the Weights and Measures Regulations. The following is a summary of the principal features only.

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SUMMARY DESCRIPTION:

CATEGORY

The device is an electronic portable bench scale (balance).

DESCRIPTION

The approved device is a self-contained electronic balance that is fitted with a single centrally located weight sensor, working on the electromagnetic force compensation principle. The weight sensor supports a circular stainless steel load receiving element. The load receiving element diameter measurements are:

<u>Model</u>	<u>Diameter</u>
AB*-S/A	84 mm
GB*-S/A	190 mm
PB*-S/A (with draft shield)	100 mm
PB*-S/A (without draft shield)	190 mm

The AB*-S/A and PB*3-S/A models have a glass draft shield as a standard feature. The other PB*-S models have the glass draft shield as an optional feature.

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The GB*-S/A and PB*-S/A models are very similar. The differences are: GB*-S/A models have no piece counting and draft shield feature. All three models: AB*-S/A, GB*-S/A and PB*-S/A are operationally identical.

INDICATOR

The device has an LCD indicator for weight display, DeltaRange for finer readability and alphanumeric messages. Annunciators are: center of zero, tare, Gross, Net, Stable, g, lb, kg.

The scale displays both “e” and “d” from no load to capacity.

FUNCTION KEYS

The device has a five button keypad that allows the following choices among others:

ON/OFF	used to turn the device on or off
O/T	used to zero the balance or to enter a tare weight
C	used to cancel functions
⇌F	used to switch between weight units.
□→	used to transfer weighing data via RS232 serial port or to calibrate or to show menu

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Note:

- Suffix -/ A indicates an approved device
- Suffix - S DR indicates "Delta Range" operation
- Suffix - S RS indicates optional RS232 communication port installed at factory

LOAD CELL CAPACITY

AB*S/A :	320 g
GB*-S/A :	8100 g
PB*-S/A :	8100 g

A mechanical overload system protects the load cells. The device can be levelled by two adjustable levelling feet and a bullseye level.

CALIBRATION

The device has an automatic calibration feature that internally re-calibrates the scale if the temperature changes enough to affect accurate weighing. This display commence to flash "autocal" when the scale senses a need to re-calibrate. Automatic calibration will only commence when the platter is empty. Calibration using external weights can be done by accessing the calibration switch.

OPERATING TEMPERATURE

AB*-S/A	12.5°C to 27.5°C
GB*-S/A	10°C to 30°C
PB*-S/A	10°C to 30°C

COMMUNICATION

S	RS232 (optional)
S	Remote printer capability (optional)

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SEALING

Means of adjustment and configuration parameters are sealed by using a wire security seal threaded through a hole in the top and bottom halves of the rear of the enclosure preventing the housing from being separated. A self-destructive seal must also be placed over one of the screws under the sub-platter and another over a hole on the rear of the scale (beneath the wall adapter input) to prevent access to the calibration switch.

EVALUATED BY

Ken Chin
Complex Approvals Examiner
Tel: (613) 952-2481
Fax: (613) 952-1754

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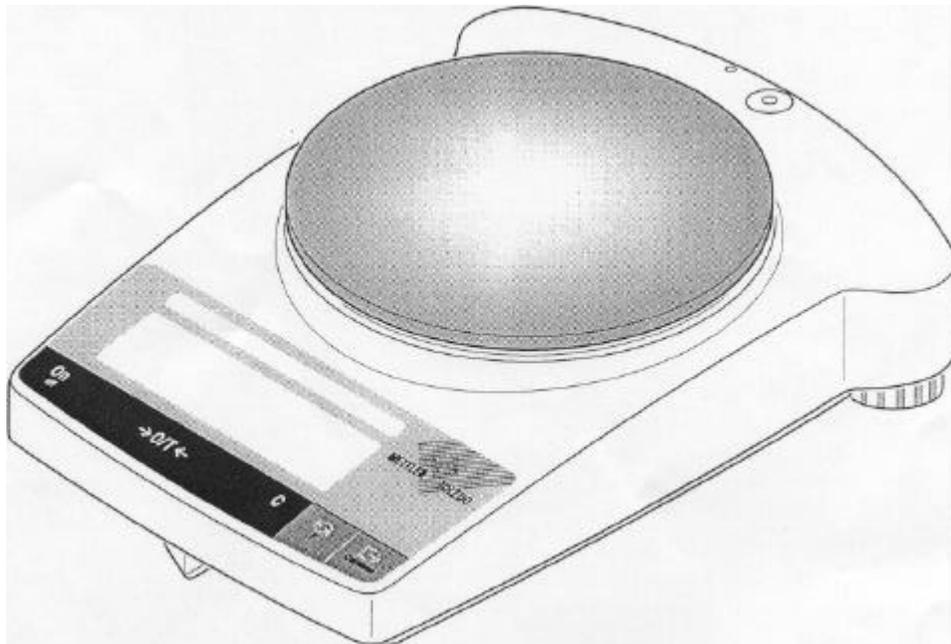
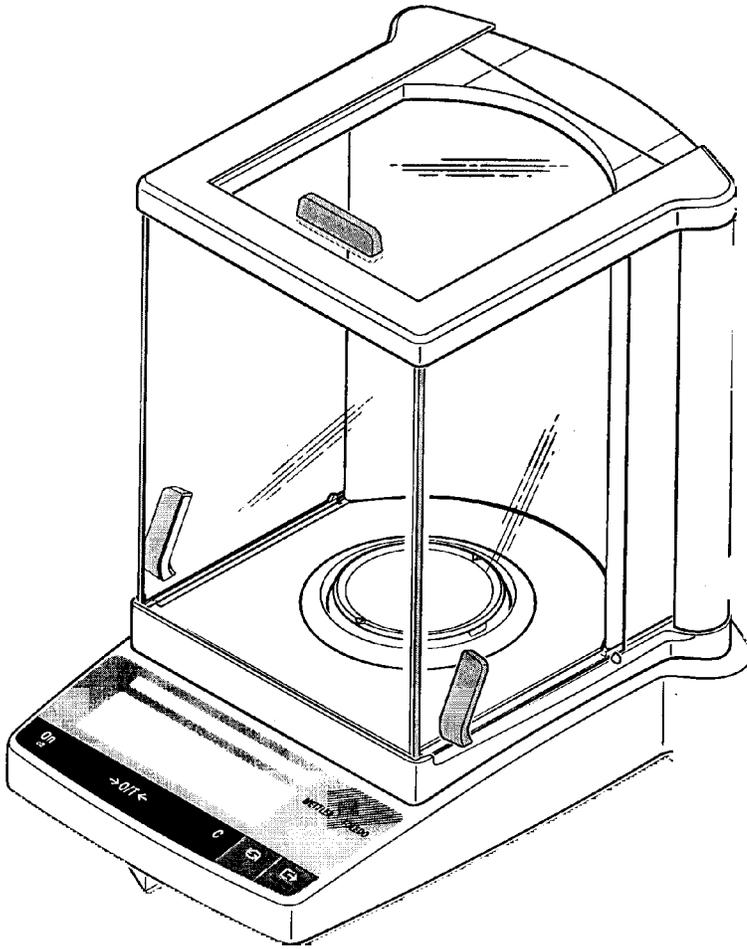
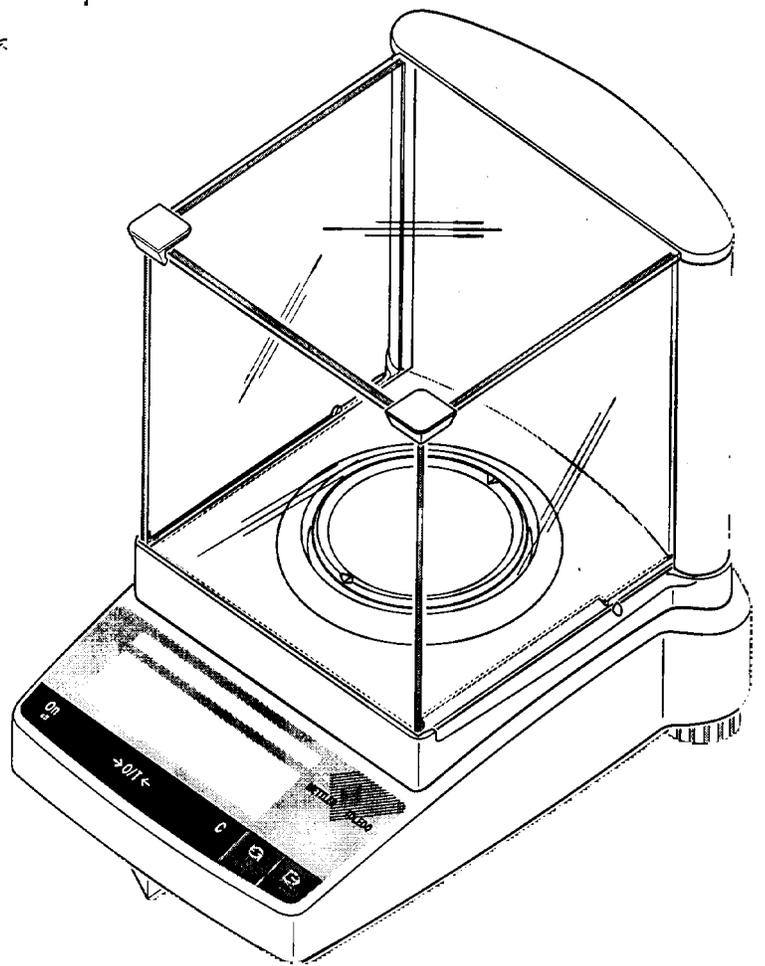
PB*-S/A

APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is hereby granted accordingly pursuant to subsection 3(1) of the said Act.

The marking, installation and manner of use of trade devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Requirements relating to marking are set forth in sections 49 to 54 of the Specifications Relating to Non-Automatic Weighing Devices. Installation and use requirements are set forth in sections 55 to 67 of the Specifications Relating to Non-Automatic Weighing Devices. A verification of conformity is required in addition to this approval. Inquiries regarding inspection and verification should be addressed to the local inspection office of Industry Canada.

Original copy signed by:

René Magnan, P. Eng
Director
Approval Services Laboratory

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